Peripex Peristaltikpumpen

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Peripex Peristaltic pumps

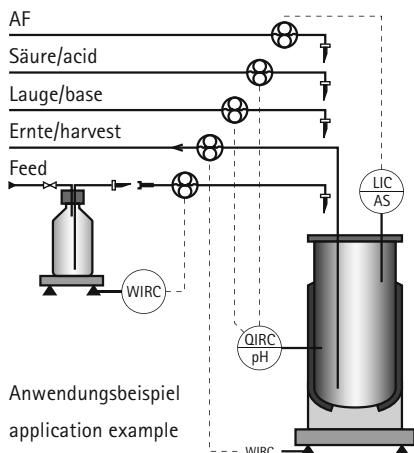
Peristaltic pumps are particularly well suitable to feed shear-sensitive, aggressive, toxic or foamy fluids. Generally they provide a good many advantages:

- easy to operate and handle
- small maintenance effort
- no contamination of the medium
- no contamination of the pump
- no dead-volume space
- self-priming.

But the most significant feature provided by the Bioengineering Peripex pumps is a precise and easy to reproduce metering and feeding. This is a consequence of the almost puls-free pumpheads line G and R and the easy to handle hoses of the pumpheads line W.

Depending on the field of application the most suitable pump can be chosen. Typical applications of Bioengineering Peripex pumps are:

- Sterile feeding and harvest in fermentation processes
- Metering of acid, base and antifoam
- Circular flow
- Continuous metered addition
- Metered addition of charges
- Sample drawing

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Type of pumps

There are two types of Bioengineering peripex pumps available:

Type 1

- Switch on/off normal pump operation, constant revolution speed
- Toggle switch to change rotational direction

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Type 2

- Integrated display and keypad
- Revolution speed is variable
- Selectable rotational direction
- Max-Speed-Function
- Profilfunktion
- Displayed consumption
- Trend displayed as graphic
- Alarm signal on display and as digital output
- Measurement and indication of consumption can be calibrated
- Control by set value and profil
- PID control of an arbitrary metered value (analog input 4...20 mA)
- Remote control of the pump by analog/digital signal (4...20 mA) or profibus
- Output of the revolution speed or the measured value with analog signal (4...20 mA) or profibus
- Settings protected by password

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Construction

The same stainless box is used for all types of pumps. A spacer plate provides the possibility to install different types of pump heads into the same box. On the rear panel the switches, connectors and interfaces are mounted. The sandwich construction makes it very easy to assembly the whole pump. The pumps can be stocked, pumps of type 2 can be arranged on edge or (as it is normal) across.

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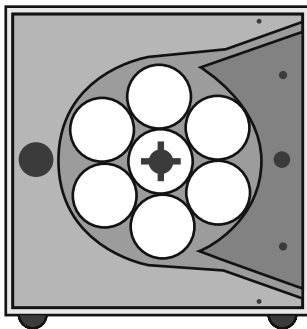
Pump heads

Line G

In this pump head's construction a new and unique principle has been applied: A central roller moves six bearing free rollers along the circular inner surface of the pump head. In this design the hose is reliably kept between the rollers and the surface. No additional clamping unit is needed. This principle and the unique form of the rollers provide a exact and practically pulsation free feeding.

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

Line R

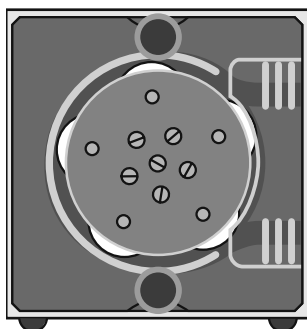This pump head is characterised by a rugged design and a very long life cycle. The twice supported mandrel drives the five rollers. The result is a precise feeding with very low pulsation. Similar to line G the mounting of the hose is very easy.

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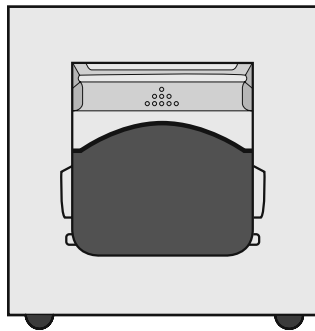The pump heads of the line G and R differ in the flow rate (see the data sheets). In a simple way they can be exchanged by each other without tools. It is possible to apply different hose dimensions.

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
Line W

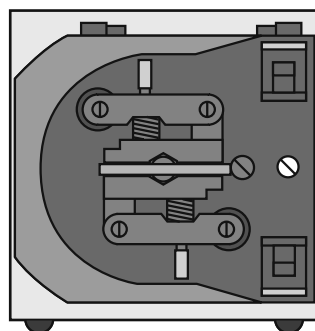A significant feature of this four roller pump head is a easy to handle device used to mount the hose in a very simple way. With a flick of the wrist the pump head can be opened and closed. An integrated device ensures the best possible clamping of the different hose sizes. In closed state the roller's area is completely protected from external influences. A bayonet socket makes it easy to disassemble the whole pump head from the casing.

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Line RL

Both spring-loaded rollers of the pumphead RL compensate automatically small deviations in the hose's wall thickness. This prolonges tube life. The handling is easy and the reproducibility high.

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For details please refer to individual data sheets

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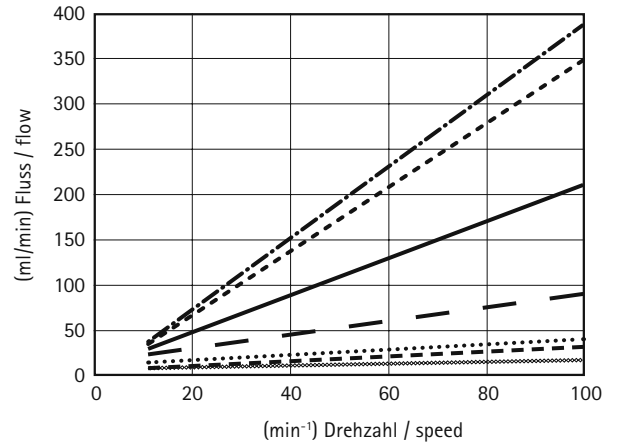
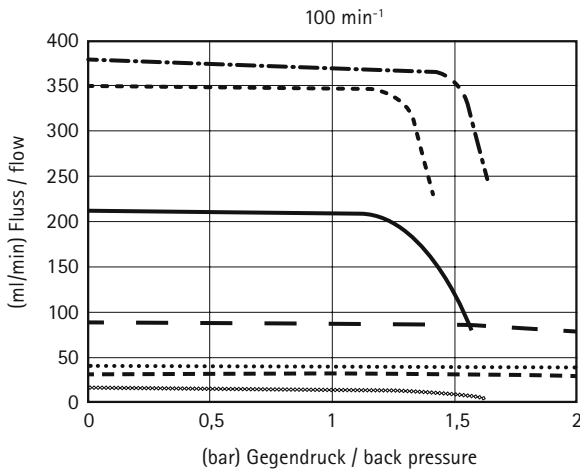
Pumpen G / R / W

Pumps G / R / W

Testmedium Wasser 20° C / Test medium water 20° C

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

Type of pump head / Type of hose

- | | |
|-----------------|--------------------|
| — — W ø 3/6 | - - - - W ø 7/10 |
| W ø 2/5,5 | - · - · - R ø 7/10 |
| - - - - G ø 3/5 | ———— W ø 5/8 |
| G ø 2/4 | |

Pumpe RL

Pump RL

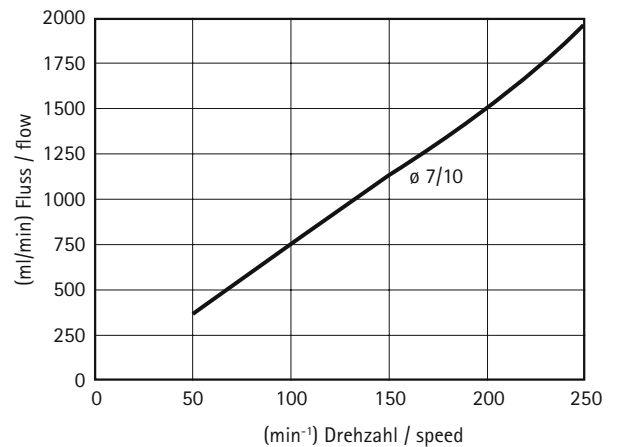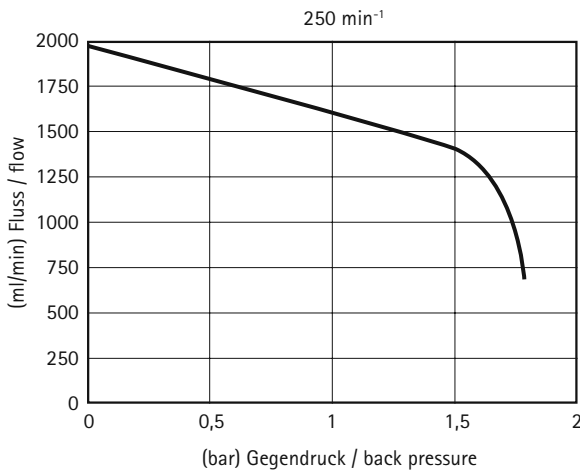Testmedium Wasser 20° C / Test medium water 20° C

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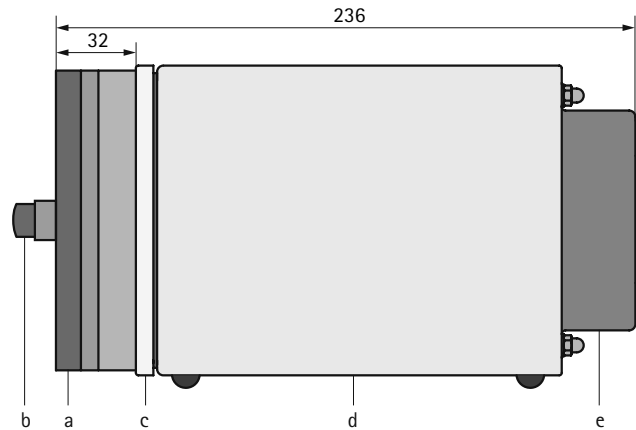
Peristaltic pump Peripex G1


- Low shear force feeding
- Very low pulsation
- Very good reproducibility
- Easy to operate
- Switch on/off
for normal pump operation
- Reversible direction of rotation

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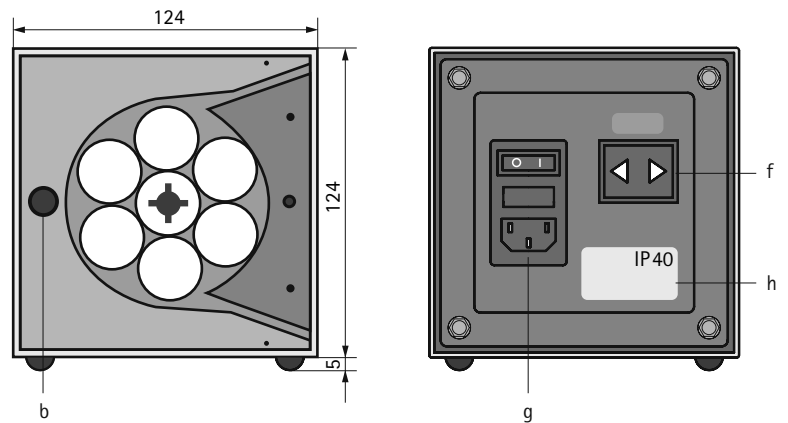
- a Pump head
- b Snap-lock to remove pump head
- c Spacer plate
- d Pump casing
- e Rear panel
- f Toggle switch for direction of rotation
- g Mains connector / power switch / fuse
- h Type plate

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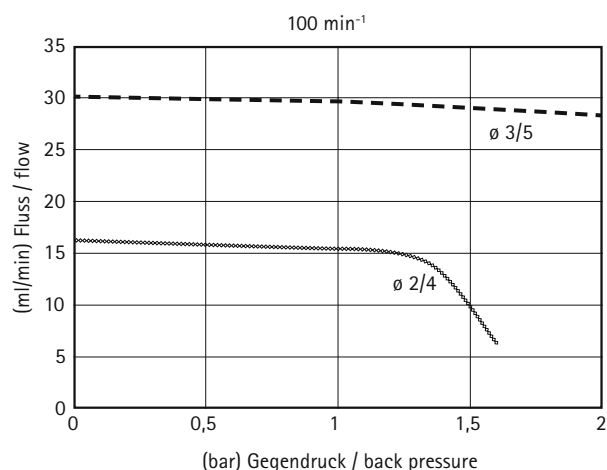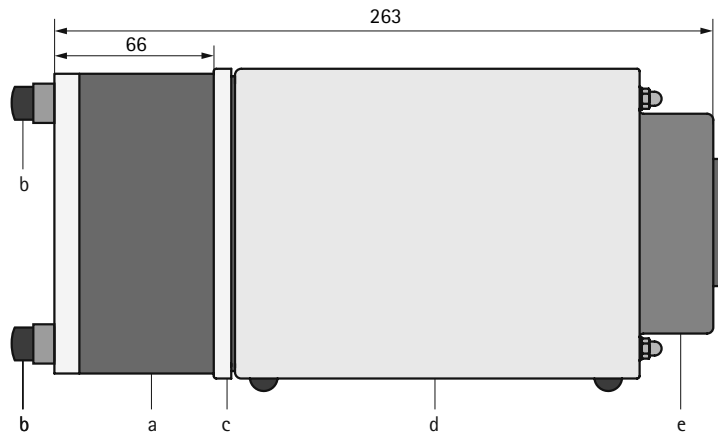
Peristaltic pump Peripex R1


- Low shear force feeding
- Very low pulsation
- Good reproducibility
- Easy to operate
- Switch on/off
for normal pump operation
- Reversible direction of rotation

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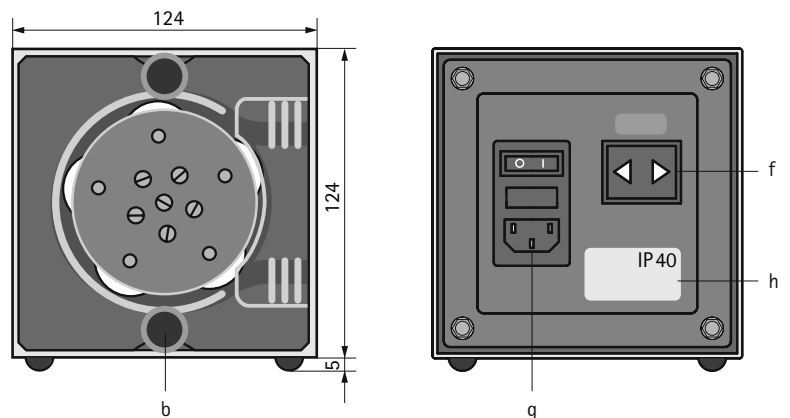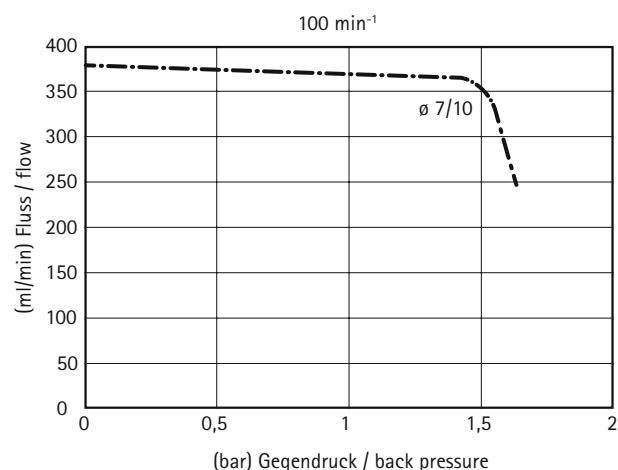
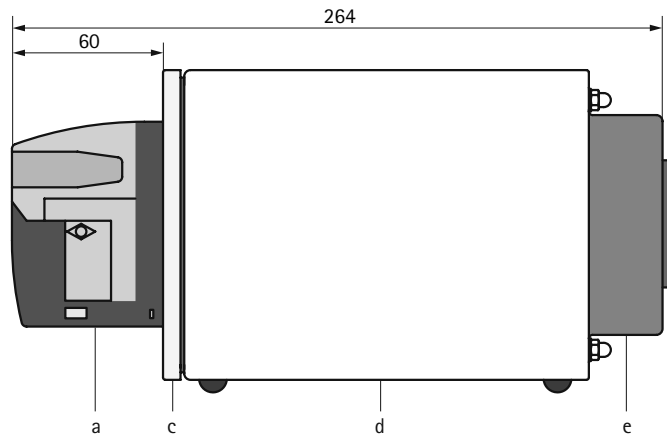
Peristaltic pump Peripex W1


- Low shear force feeding
- Low pulsation
- Good reproducibility
- Easy to operate
- Switch on/off
for normal pump operation
- Reversible direction of rotation

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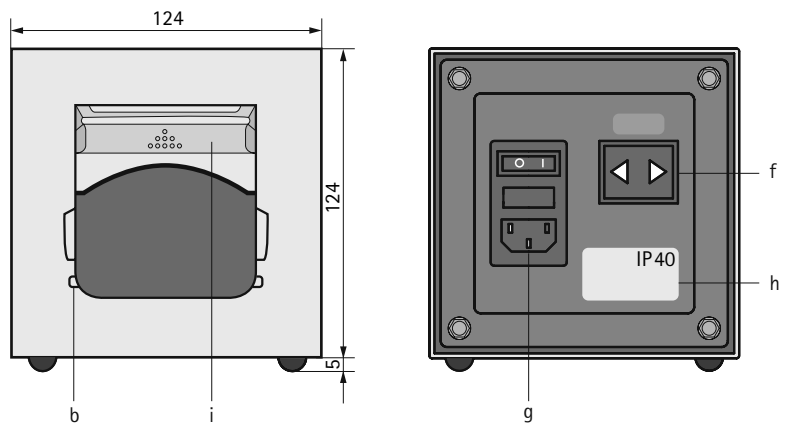
- a Pump head
- b Hose clamp
- c Spacer plate
- d Pump casing
- e Rear panel
- f Toggle switch for direction of rotation
- g Mains connector / power switch / fuse
- h Type plate
- i Cap

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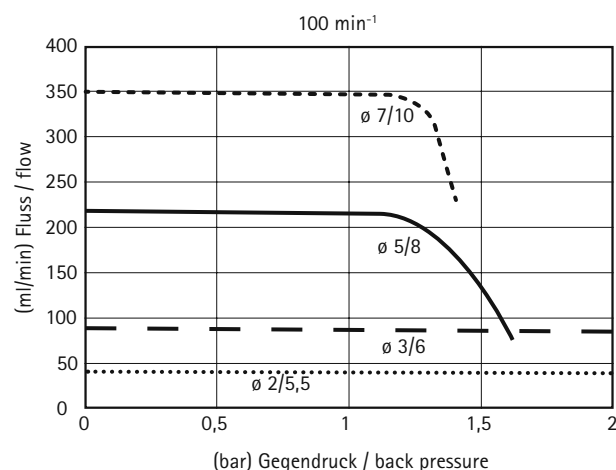
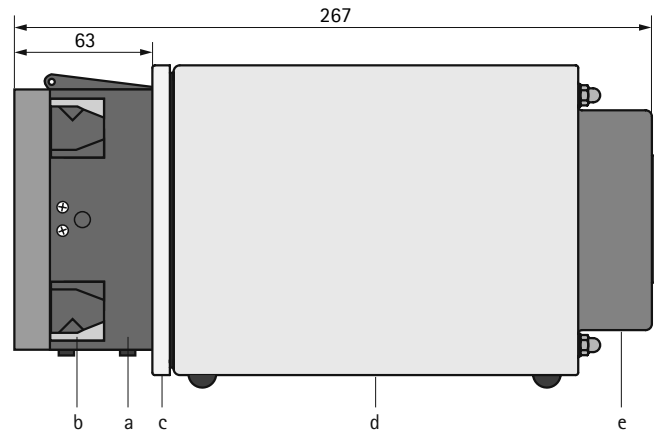
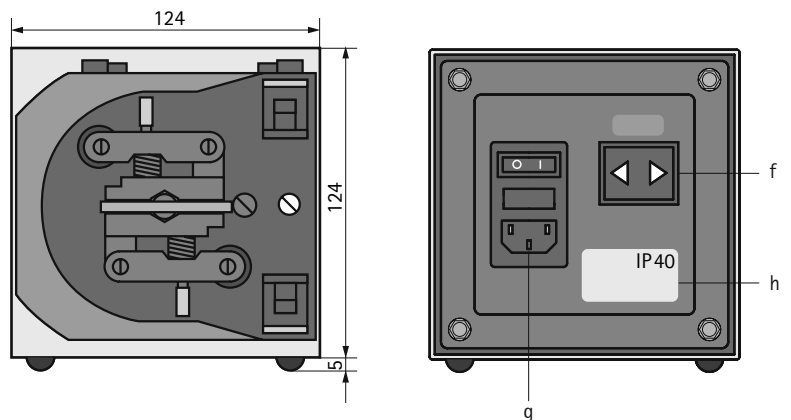
- a Pump head
- b Hose clamp
- c Spacer plate
- d Pump casing
- e Rear panel
- f Toggle switch for direction of rotation
- g Mains connector / power switch / fuse
- h Type plate

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
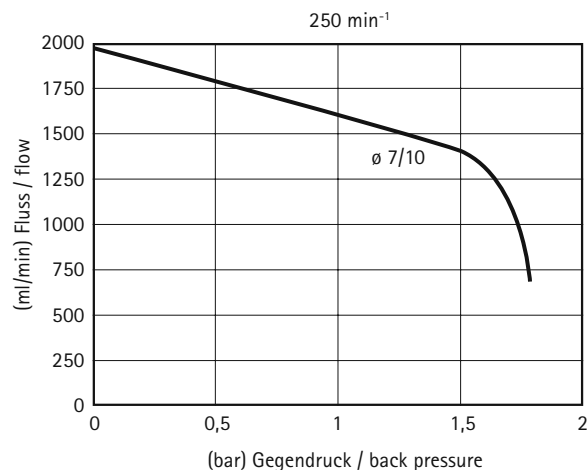
Peristaltic pump Peripex G2 IP40

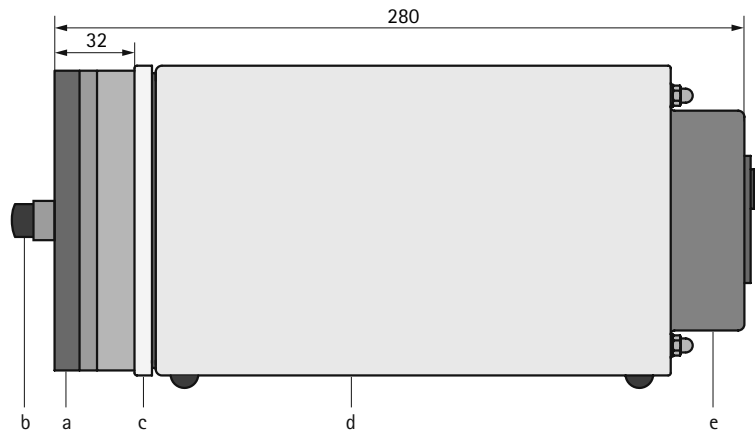
- Low shear force feeding
- Very low pulsation
- Very good reproducibility
- Operated by keypad functions
- Display for status indication and parameter input
- Reversible direction of rotation

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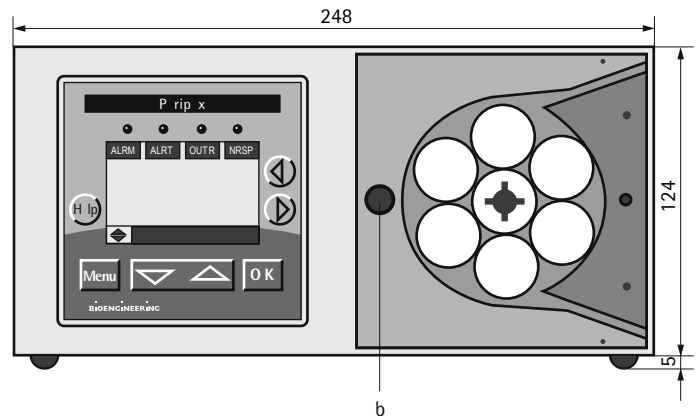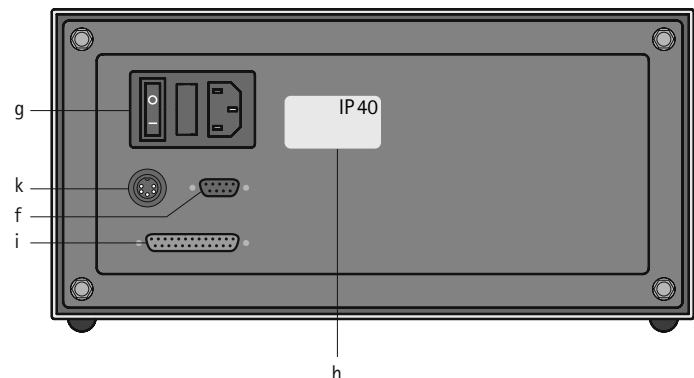- a Pump head
- b Snap-lock to remove pump head
- c Spacer plate
- d Pump casing
- e Rear panel
- f Profibus DP
- g Mains connector / power switch / fuse
- h Type plate
- i Remote control connector
- k Cascade

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

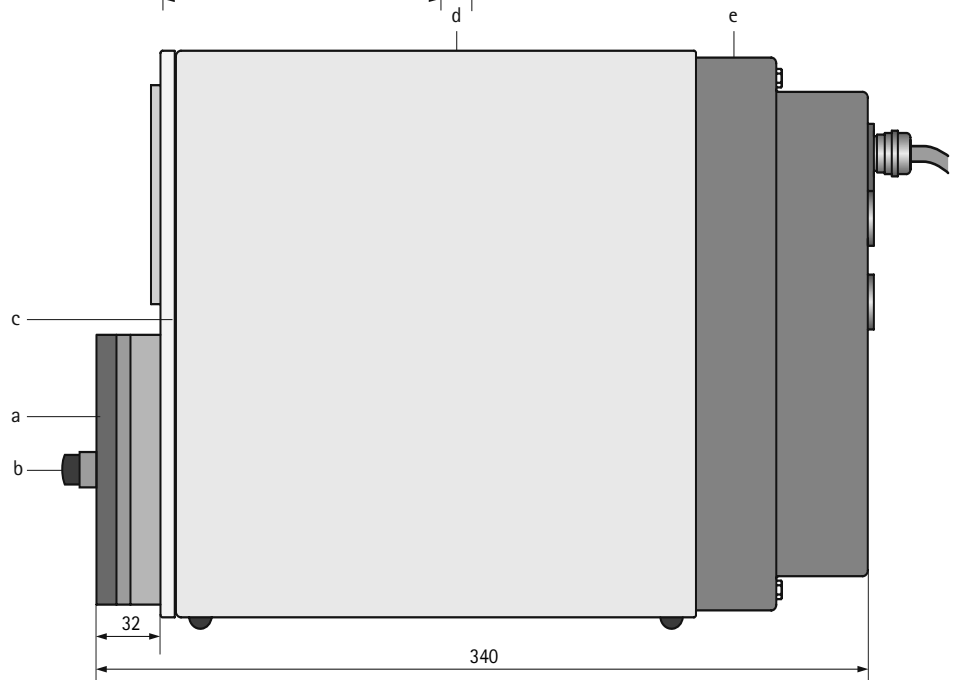
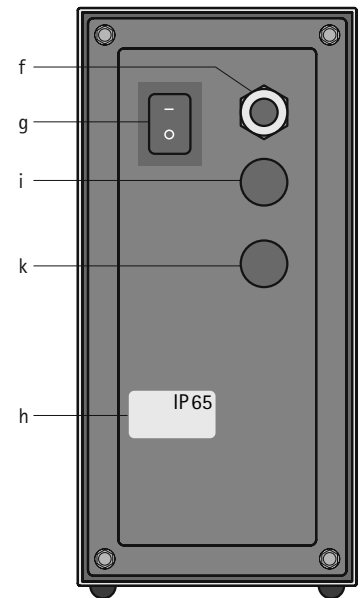
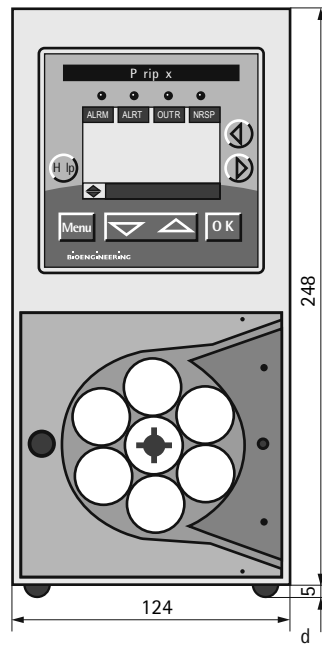
- a Pump head
- b Snap-lock to remove pump head
- c Spacer plate
- d Pump casing
- e Rear panel
- f Mains connector
- g Power switch
- h Type plate
- i For cable clamp
- k For cable clamp

Technische Daten G2	Technical data G2	
(Technische Änderungen vorbehalten)	(Subject to technical modifications)	
Pumpenkopf	Pump head	G
Anzahl Rollen	Number of rolls	6
Min. Schlauchinnen \varnothing	Hose inner \varnothing , min.	1,15 mm
Max. Schlauchinnen \varnothing	Hose inner \varnothing , max.	3 mm
Schlauchwandstärke	Hose wall thickness	1 mm
Drehzahl	Revolutions	0-100 min ⁻¹
Reproduzierbarkeit ohne Gegendruck	Reproducibility without back pressure	$\pm 1\%$
Motortyp DC Getriebemotor	Gear motor	24V, DC
Leistungsaufnahme	Power consumption	25 W
Spannung	Line voltage	100...230V, AC
Frequenz	Line frequency	50...60 Hz
Schutzart	Protection class	Standard IP40
Schutzart optional	Protection class, optional	IP65
Gewicht IP40 / IP65	Weight IP40 / IP65	6,4 kg / 6,7 kg
Gehäusematerial	Housing material	Inox, geschliffen Inox, grinded
Rückblech	Rear panel	Stahl, pulverbeschichtet RAL 5010 Steel, powder coating RAL 5010
Pumpenkopfmateri	Material of pumphead	Acrylglas, polymethacrylate

Bestellangaben	Ordering information	Art. Nr. / Part no.
Peripex G2 horizontal IP40 (Standard)	Peripex G2 horizontal IP40 (Standard)	13935.1
Peripex G2 vertikal IP40	Peripex G2 vertical IP40	13935.3
Peripex G2 horizontal IP65	Peripex G2 horizontal IP65	13934.1
Peripex G2 vertikal IP65	Peripex G2 vertical IP65	13934.3
Schlauch $\varnothing 2/4$ Silikon transparent	Hose $\varnothing 2/4$ silicone transparent	53444
Schlauch $\varnothing 3/5$ Silikon transparent	Hose $\varnothing 3/5$ silicone transparent	50222
Pumpenkopf G	Pump head G	53084
kompatibel zu Peripex R	compatible with Peripex R	
Zubehör	Accessories	Netzkabel IP40 / Power cord IP40

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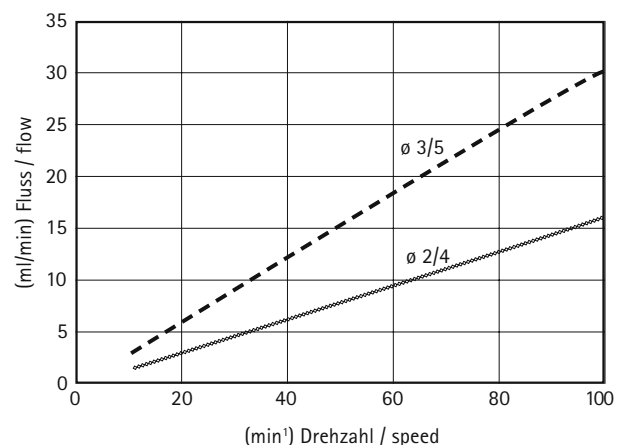
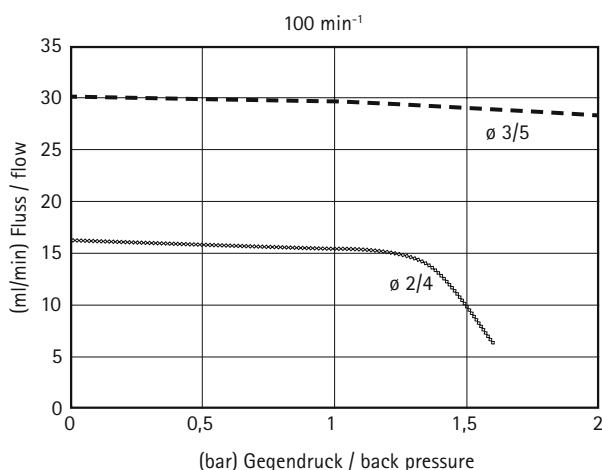
Peristaltic pump Peripex R2 IP40

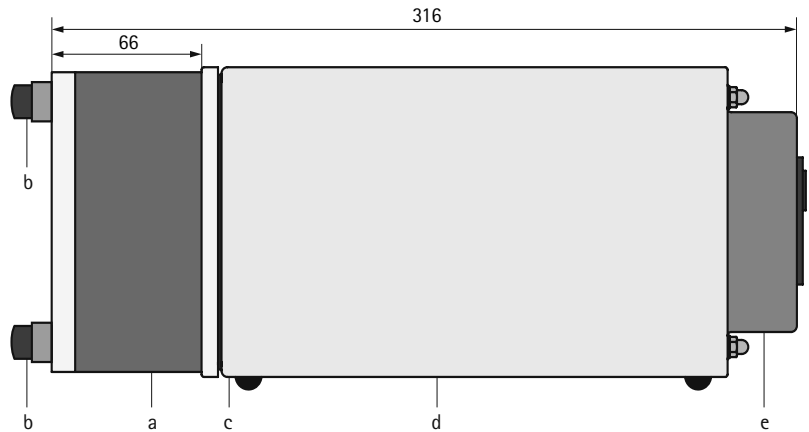
- Low shear force feeding
- Very low pulsation
- Good reproducibility
- Operated by keypad functions
- Display for status indication and parameter input
- Reversible direction of rotation

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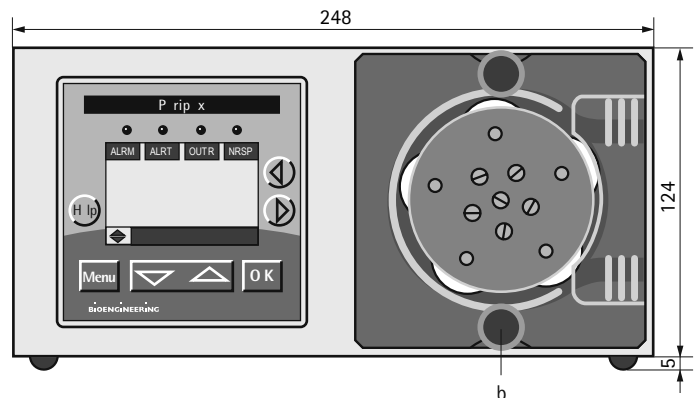
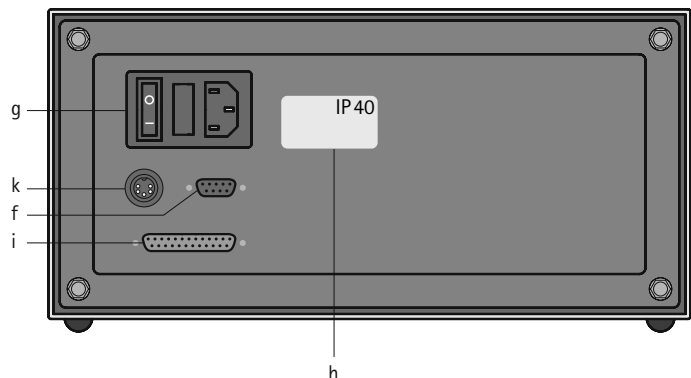
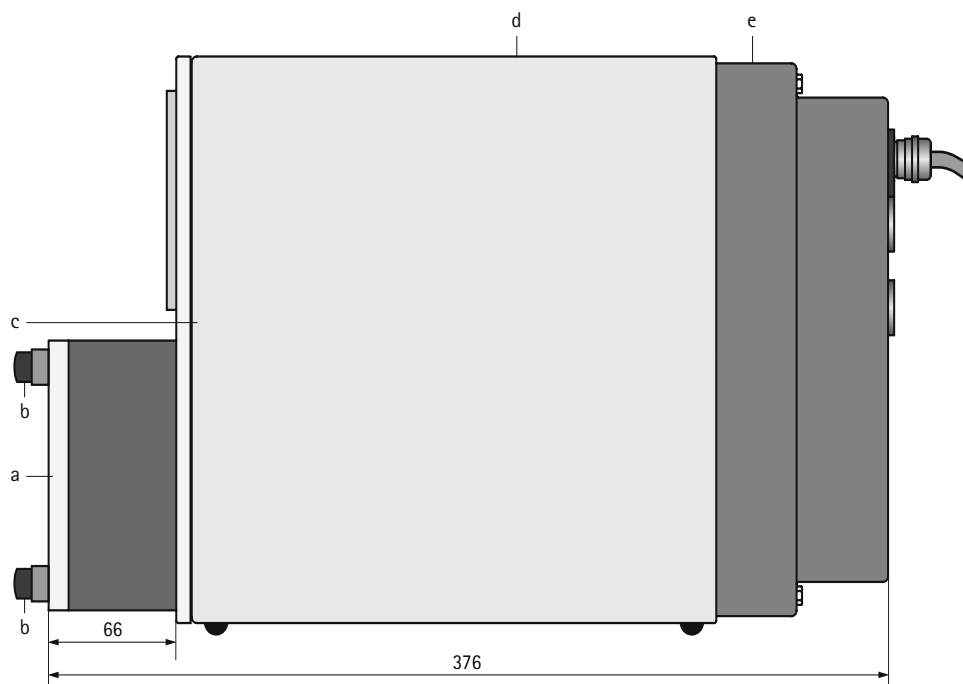
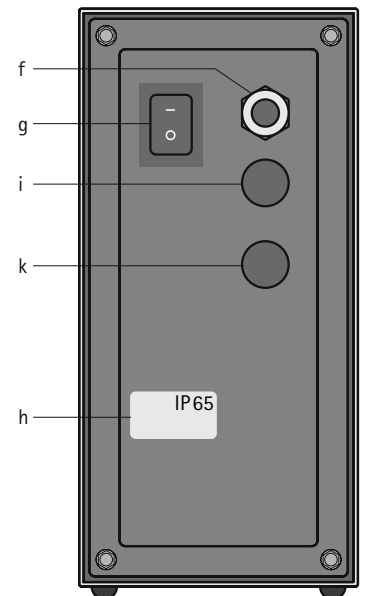
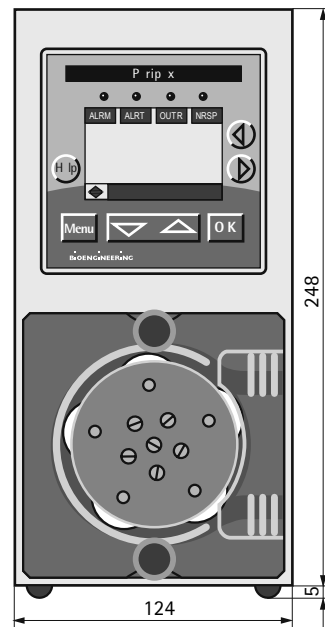
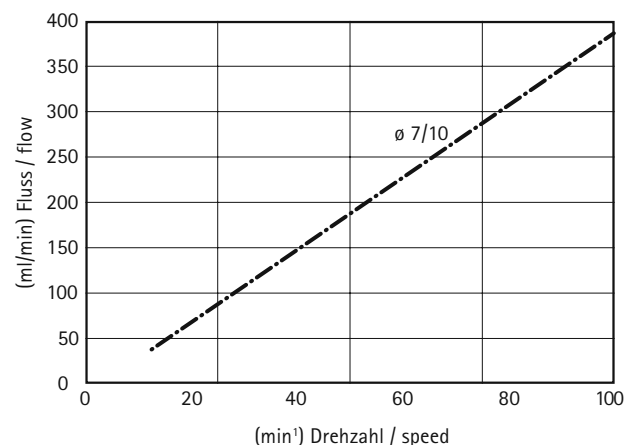
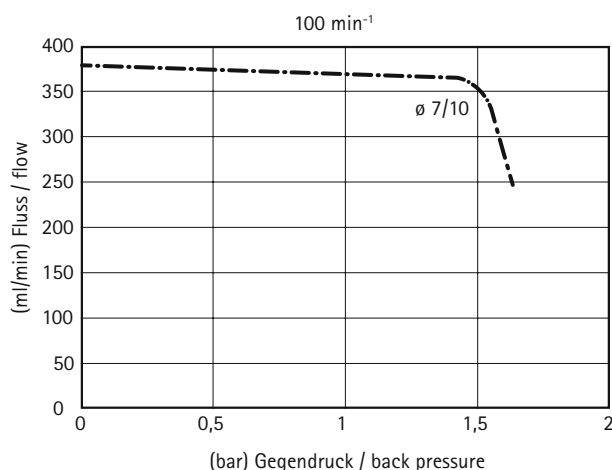
Peristaltic pump Peripex W2 IP40

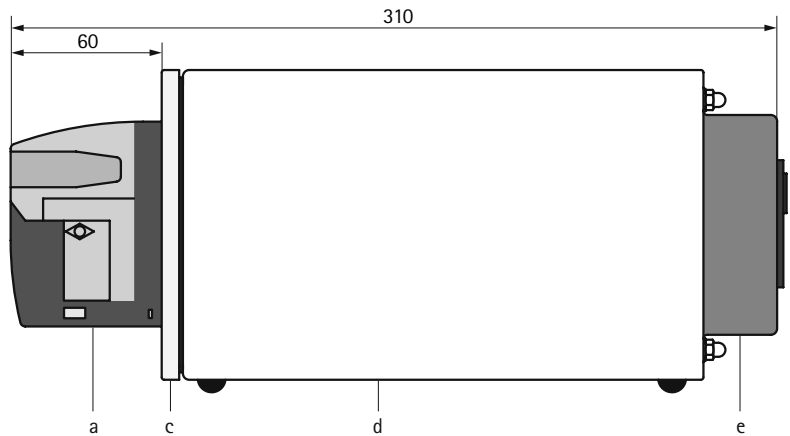
- Low shear force feeding
- Low pulsation
- Good reproducibility
- Operated by keypad functions
- Display for status indication and parameter input
- Reversible direction of rotation

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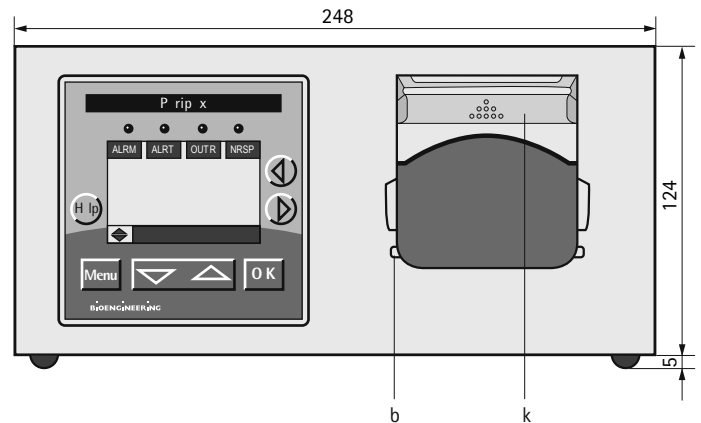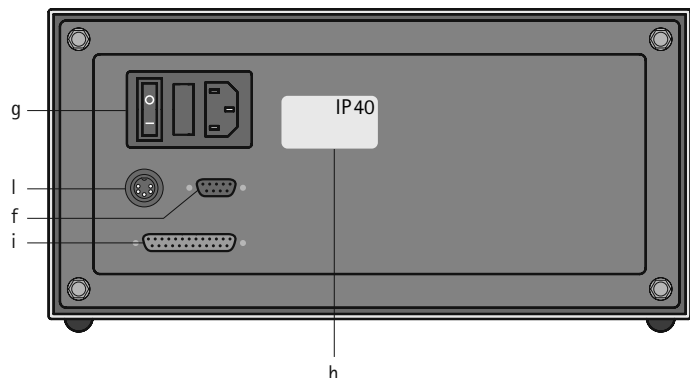- a Pump head
- b Hose clamp
- c Spacer plate
- d Pump casing
- e Rear panel
- f Profibus DP
- g Mains connector / power switch / fuse
- h Type plate
- i Remote control connector
- k Cap
- l Cascade

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
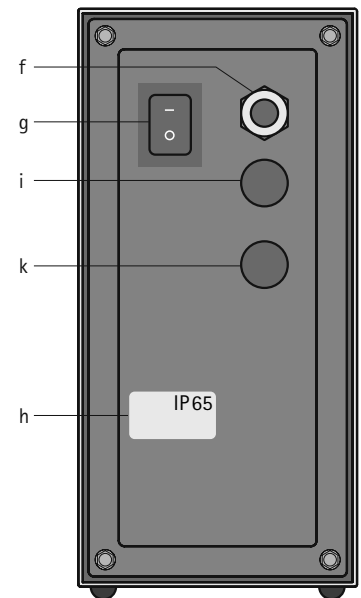
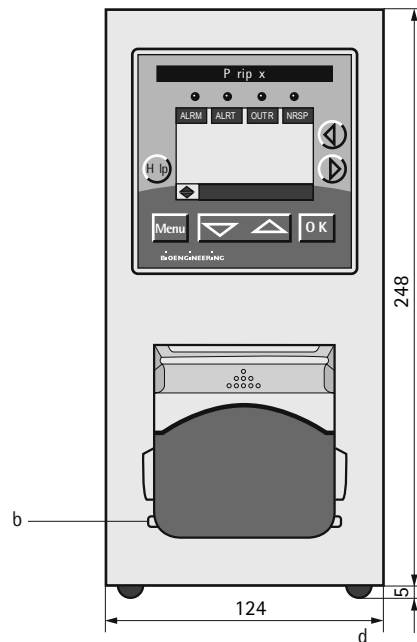
- Low shear force feeding
- Low pulsation
- Good reproducibility
- Operated by keypadfunctions
- Display for status indication and parameter input
- Reversible direction of rotation

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
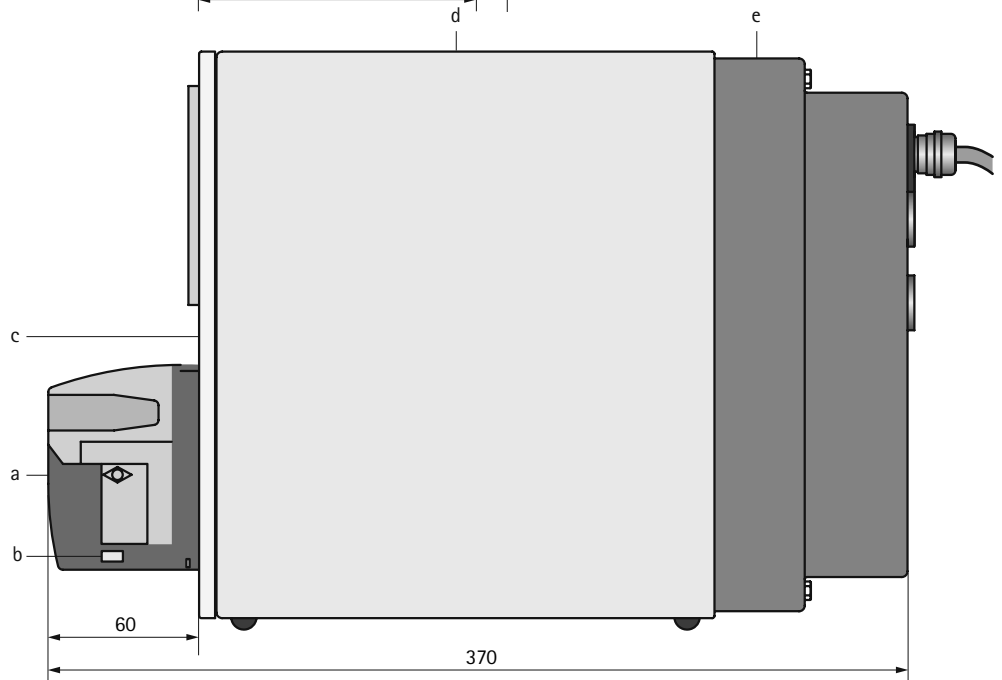
- a Pump head
- b Hose clamp
- c Spacer plate
- d Pump casing
- e Rear panel
- f Mains connector
- g Power switch
- h Type plate
- i For cable clamp
- k For cable clamp

Technische Daten W2	Technical data W2	
(Technische Änderungen vorbehalten)	(Subject to technical modifications)	
Pumpenkopf	Pump head	W
Anzahl Rollen	Number of rolls	4
Min. Schlauchinnen \varnothing	Hose inner \varnothing , min.	0,5 mm
Max. Schlauchinnen \varnothing	Hose inner \varnothing , max.	8 mm
Schlauchwandstärke	Hose wall thickness	1,6 mm
Drehzahl	Revolutions	0-100 min ⁻¹ var
Reproduzierbarkeit ohne Gegendruck	Reproducibility without back pressure	
Schlauch \varnothing 2/5,5	Hose \varnothing 2/5,5	\pm 1 %
Schlauch \varnothing 3/6	Hose \varnothing 3/6	\pm 1 %
Schlauch \varnothing 5/8	Hose \varnothing 5/8	\pm 0,5 %
Schlauch \varnothing 7/10	Hose \varnothing 7/10	\pm 0,5 %
Motortyp DC Getriebemotor	Gear motor	24V, DC
Leistungsaufnahme	Power consumption	25 W
Spannung	Line voltage	100...230V, AC
Frequenz	Line frequency	50...60 Hz
Schutzart	Protection class	Standard IP40
Schutzart optional	Protection class, optional	IP65
Gewicht IP40 / IP65	Weight IP40 / IP65	6,1 kg / 6,4 kg
Gehäusematerial	Housing material	Inox, geschliffen / grinded
Rückblech		Stahl, pulverbeschichtet RAL 5010
	Rear panel	Steel, powder coating RAL 5010
Pumpenkopfmaterial	Material of pumphead	Acrylglas, polymethacrylate

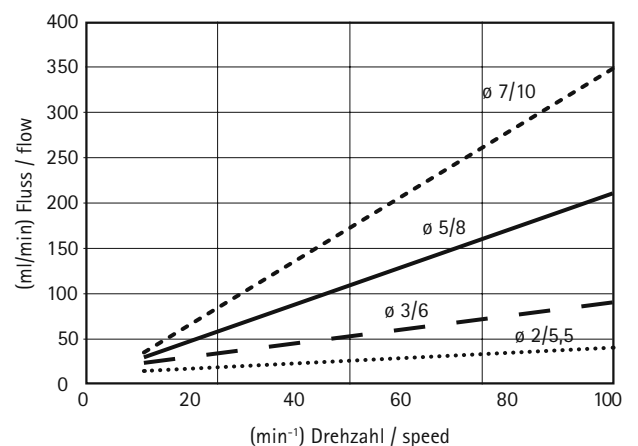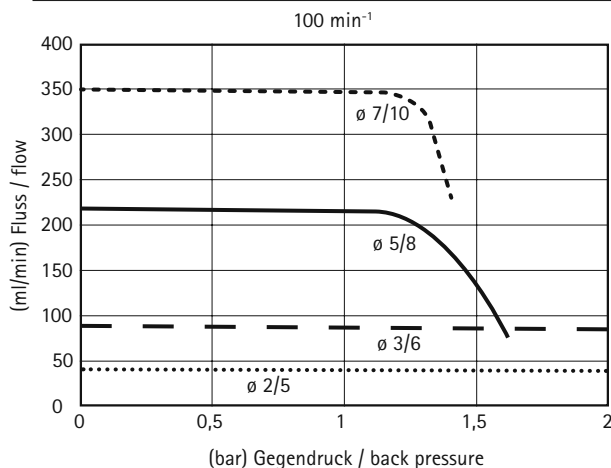Bestellangaben	Ordering information	Art. Nr. / Part no.
Peripex W2 horizontal IP40 (Standard)	Peripex W2 horizontal IP40 (Standard)	13936.1
Peripex W2 vertikal IP40	Peripex W2 vertical IP40	13936.2
Peripex W2 horizontal IP65	Peripex W2 horizontal IP65	13937.1
Peripex W2 vertikal IP65	Peripex W2 vertical IP65	13937.2
Schlauch \varnothing 2/5,5 Silikon transparent	Hose \varnothing 2/5,5 silicone transparent	51808
Schlauch \varnothing 3/6 Silikon transparent	Hose \varnothing 3/6 silicone transparent	53446
Schlauch \varnothing 5/8 Silikon transparent	Hose \varnothing 5/8 silicone transparent	53447
Schlauch \varnothing 7/10 Silikon transparent	Hose \varnothing 7/10 silicone transparent	50224
Pumpenkopf W	Pump head W	53443
Zubehör	Accessories	Netz Kabel IP40 / Power cord IP40

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